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An important property of liquids is their viscosity. A measure for the viscosity is the flow time, determined at a specified temperature, using a flow cup.

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## Application areas

- for paint and varnish manufacturers, its customers and test laboratories
- the test can be carried out in the lab and during production

## Standard delivery

- 1 flow cup/immersion flow cup
- 1 certificate of manufacturer
- 1 storage box

## Options

- ZTJ 3020 Temperature control jacket
- ACC050 insertion thermometer up to 400°C (752°F) (for fluids)
- ACC072 three-foot stand completely made of stainless steel, levelling, adjustable feet, with spirit level, height of 260 mm (10.24")
- ACC152 glass plate
- ZKZ005 calibration and certification (incl. certificate)

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## Handling flow cups

- choose a flow cup that will give a flow time of between 30 s and 100 s (depending on the standard)
- put the flow cup into the levelled three-foot stand
- close the orifice of the cup by a finger and fill the cup with the liquid under test, avoiding the formation of air bubbles
- slide a glass plate over the rim of the cup so that no air bubbles will be formed between the glass plate and the surface of the sample
- draw the glass plate horizontally away, remove the finger from the orifice of the cup and simultaneously start the timing-device
- stop the timing-device as soon as the stream of the sample first breaks close to the orifice

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## Handling immersion flow cups

- immerse the immersion flow cup entirely below the surface of the liquid under test
- withdraw the cup vertically and start the timing-device
- stop the timing-device as soon as the stream of the sample first breaks close to the orifice
- the average out of 3 measurements will give the flow time

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## Technical specification

version	cup	with fixed jet*	standard	image
ZFC 3010	flow cup (Ford)	No. 1, 2, 3, 4 or 5.	ASTM D 1200	
ZFC 3011	flow cup	No. 1, 2, 3, 4, 5, 6, 7 or 8.	**	
ZFC 3012	immersion flow cup	No. 1, 2, 3, 4, 5, 6, 7 or 8.	**	
ZFC 3013	flow cup	No. 3, 4, 5 or 6	ISO 2431	
ZFC 3014	immersion flow cup	No. 3, 4, 5 or 6	ISO 2431	
ZFC 3015	immersion flow cup (Ford)	No. 1, 2, 3, 4 or 5.	ASTM D 1200	

\* Please specify jet No. when ordering.  
 \*\* Jet No. 4 in acc. with DIN 53211, withdrawn in October 1996  
 material: cup in anodised aluminium, jet in stainless steel  
 warranty: 2 years