Metallpyknometer

Metal pycnometer

ZPM 3030

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The density is an essential quality characteristic, for example it allows recognising errors regarding the composition and gives indications of the purity of a product.

The test principle is based on the weighing of a given volume of the liquid under test, as it is defined by the pycnometer. From the difference of the masses of the filled and the empty pycnometer as well as the known filling volume the density is calculated.

Application areas

- for very different branches of industry, e.g. for the paint, cosmetics, food, cleaning agents industries and the chemical industry in general, as well as for technical high schools, universities and testing institutes
- laboratory apparatus for the quality control and for research and development

Standard delivery

- 1 pycnometer, consisting of a cylindrical beaker with a lid having a hole
- 1 certificate of manufacturer
- 1 certificate of calibration
- 1 storage box

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Handling

- weigh the carefully cleaned empty pycnometer plus lid and record the mass
- bring the pycnometer and the substance under test to the temperature required in the chosen standard unless otherwise agreed
- fill the pycnometer with the substance under test and place the lid firmly in position without tilting, whilst avoiding inclusion of air bubbles
- remove any excess liquid from the lid using a clean and absorbent cloth
- weigh the filled pycnometer
- calculate the density, using the following formula:
density = mass (g) of filled pycnometer minus mass (g) of empty pycnometer, divided by the volume (ml) of the pycnometer
- do not use sharp-edged objects when cleaning the pycnometer in order to avoid damaging, but a suitable solvent

Technical specification

version	filling volume	height/diameter	weight
ZPM 3030.50	50 ml	35 mm x Ø 54 mm (1.38" x Ø 2.13")	144.1 g (0.318 lbs)
ZPM 3030.100	100 ml	63 mm x Ø 54 mm (2.48" x Ø 2.13")	174.4 g (0.384 lbs)

material: stainless steel
 manufacturing tolerance: ± 0,2 %, related to the volume; as option with certificate of calibration ± 0,1 %
 standards: ASTM D 1475, EN ISO 2811-1, SNV 37100, VDA 621-103
 warranty: 2 years